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(a) Species Specific PCR (mitochondrial D Loop)

(b) PCR by Universal Primer Technique (12s rRNA) & Species Identification by RFLP

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Standard Operative Procedure for Sampling, Handling, Handover & Documentation of Bovine Tissues collected for analysis in Forensic Analysis Laboratory for Detection of Bovine Tissue :

To ensure protection of cow progeny, prohibit slaughter & smuggling of cow progeny, Government Uttarakhand has enacted '**Uttarakhand Protection of Cow Progeny, Act 2007**'. Under '**section 13(1)**' of the act, '**Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018**' have been framed.

Under '**Rule-33**' of the '**Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018**', Government Uttarakhand has issued notification No.-1002/ XV-1/ 19/7(14)/2005 Pashupalan Section-1, dated 1st November, 2019 for, establishment of 'Forensic Analysis Laboratory for Detection of Bovine Tissue'. 'Forensic Analysis Laboratory for Detection of Bovine Tissue' has now been established at Sheep & Wool Extension Institute, Pashulok Rishikesh, Dehradun.

In the light of '**Rule-34**' under the '**Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018**' the Standard Operating Procedure for collection, sealing, handling, transport, receiving, review, acceptance/ rejection, analysis & declaration of results of analysis of the suspected Bovine Tissue Sample is being prescribed as following :-

1. Project Director, Sheep & Wool Extension Institute, Pashulok Rishikesh, Dehradun, shall be Head/ In-charge of the Bovine Tissue Forensic Analysis Laboratory, having Financial and Administrative control of the Laboratory. Project Director Pashulok shall be doing all the communication & release of Laboratory Analysis Report on behalf of the 'Forensic Analysis Laboratory for Detection of Bovine Tissue'.
2. On receiving the request on '**PROFORMA-17 (under Rule-34 of Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018)**', from Controlling Officer of the Police Thana/ Police Chauki of the area, Jurisdictional Veterinary Officer of Government Veterinary Hospital of the area shall collect & preserve the Suspected Bovine Tissue Samples, get it signed by Police Personnel & self, seal the Sample and hand over to the Police Personnel, along with the sample seal on a paper inside envelop and complete documentation as prescribed in '**PROFORMA-18 under Rule-34 of Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018**'.

3. Jurisdictional Veterinary Officer of Government Veterinary Hospital of the area shall ensure following precautions while collecting & preserving the Suspected Bovine Tissue Samples :-
- A. Preferably uncooked flesh sample should be collected, if not available, then different type of organs or tissues may be collected. Better to avoid the sampling of putrefied tissues. DNA may not get recovered from putrefied sample, barring the identification of bovine tissue while laboratory analysis.
 - B. 100g of sample should be forwarded to Project Director, Pashulok. Samples in insufficient quantity shall not be accepted.
 - C. Before sealing, sample should be collected in Sample Container Box with proper label narrating the type of tissue.
 - D. Sample should be preserved & completely covered from all the sides, in the COMMON SALT while putting it in Sample Container Box. Thread used to pack wrap the Sample Container Box should be properly sealed with lacquer seal. Sample Container Box should be packed in cloth and the sewing on the cloth should be properly sealed. Preferably Sample Container Box should be maintained in cold chain, so that it may not get putrefied.
 - E. Samples **preserved or contaminated with FORMALIN shall not be** accepted by Project Director Pashulok.
 - F. Flesh/ Tissue identified in different containers/ pouches should be collected in different sample containers with different labels.
 - G. Sample Tissue collected in one sample container shall be considered as one sample.
 - H. **Samples without intact lacquer seal shall not be accepted by Project Director Pashulok.**
 - I. Jurisdictional Veterinary Officer shall collect & seal the Sample in presence of representative from police department. After getting the slips signed & sealed by representative from police department and self, the documentation shall be done as per the PROFORMA-18 (*under Rule-34 of Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018*). Sealed Sample, Sample Seal put on a paper inside envelop and complete documentation as prescribed in **PROFORMA-18** shall handed over to representative from police department. After completion of documentation at the level of Police

Thana/ Police Chauki, the same may be promptly delivered to Project Director (Ex-officio In-charge, Forensic Analysis Laboratory for Detection of Bovine Tissue), Pashulok Rishikesh, Dehradun.

- J. Jurisdictional Veterinary Officer shall attest the Sample Seal put on a Paper, put it in envelope & handover to Representative from Police Department.
4. After receiving the sample, Controlling Officer of the Police Thana/ Police Chauki of the area, shall complete the entry of details in **PROFORMA-18** (prescribed as per '**Rule-34**' under the '**Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018**') & promptly dispatch the sealed samples under the police custody along with the following documents :-
 - A. PROFORMA-18 signed by sample submitting authority.
 - B. Copy of the FIR (incase FIR been filed) / brief detail of complaint.
 - C. Sample should properly sealed with lacquer along with the signature slips signed by Jurisdictional Veterinary Officer & Representative from Police Department.
 - D. Attested copy of the Departmental Identity Card of the police personnel accompanying the sample.
 - E. Envelope containing paper with Sample Seal attested by Jurisdictional Veterinary Officer.
 5. Project Director Pashulok shall be accepting the Samples only under the condition that, the above mentioned norms prescribed in Point No. 3 & 4 are properly fulfilled.
 6. Samples dispatched by courier/ postal mail shall not be accepted by Project Director Pashulok.
 7. Sample submitted to Project Director Pashulok, shall be opened and reviewed in presence of sample accompanying police personnel and Head/ In-charge of the Laboratory, as prescribed in 'Format-1'.
 8. Samples found acceptable while review, shall be issued the Receipt by Project Director Pashulok, as prescribed in 'Format-2'. In case samples found not fit to be accepted for further analysis, shall be rejected and all the concerned, shall be informed on the format as prescribed in 'Format-3'. Copy the same shall be shared to concerning Veterinary Officer (who sealed the sample), Chief Veterinary Officer of the concerning district & undersigned

9. Putrefied samples may not get detected while laboratory analysis because of degenerated DNA.
10. After accepting the sample, Project Director Pashulok shall ensure entry of details in **Sample Receiving Register** as prescribed in 'Format-4'. Each Sample shall be given an Identity Number. This Identity Number shall be used for further Laboratory Analysis/ Storage of the samples. Only after labeling of each Sample with it's Identity Number, the sample shall be handed-over for further analysis in the Forensic Laboratory.
11. Project Director Pashulok shall ensure the confidentiality about origin of samples with the Laboratory officials/ helping staff involved in Laboratory Analysis/ Storage of the samples. Project Director Pashulok shall keep the **Sample Receiving Register** in his custody.
12. As prescribed on 'Format-5', Project Director Pashulok shall hand-over each of the Sample with Individual Identity Number, to Veterinary Officer Grade-1/ Veterinary Officer Grade-2 of the Laboratory for further analysis.
13. Veterinary Officer Grade-1/ Veterinary Officer Grade-2 of the Laboratory shall preserve & seal the half quantity of each of the sample for storage up to one year & use remaining half of the sample for lab analysis.
14. Standard Operative Procedure for following two techniques has been approved for laboratory analysis of the samples :-
 - (a) Species Specific PCR (mitochondrial D Loop)
 - (b) PCR by Universal Primer Technique (12s rRNA) & Species Identification by RFLP
15. Veterinary Officer Grade-1/ Veterinary Officer Grade-2 of the Laboratory shall do analysis of the samples as per the Standard Operative Procedure approved by undersigned.
16. After analysis of the samples, Veterinary Officer Grade-1/ Veterinary Officer Grade-2 shall enter the report in **Sample Analysis Register of the Laboratory** as prescribed in 'Format-6' and submit the same to Project Director Pashulok Rishikesh (Head/ In-charge of the Bovine Tissue Forensic Analysis Laboratory).
17. Project Director Pashulok Rishikesh shall convey the signed & sealed copy of the **Sample Analysis Report** on 'Format-7' to the Controlling Officer of the concerning Police Thana/ Police Chauki. Copy of the same Report shall be shared to concerning Veterinary Officer (who sealed the sample), Chief Veterinary Officer of the concerning district & undersigned.
18. In case needed, Bovine Tissue Forensic Analysis Laboratory may get time to time guidance/ help/ assistance/ skill improvement/ validation from higher

& better equipped laboratory of the Research Institutions functioning under Government of India.

19. To ensure better quality control, fairness & revalidation of operations, randomly remaining part of the Stored Sample may be sent to higher & better equipped laboratory of the Research Institutions functioning under Government of India for reconfirmation of the results.
20. The Sample Analysis Report will be submitted to the forwarding authority within two months.

**FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE
SHEEP AND WOOL EXTENSION INSTITUTE
Pashulok, Rishikesh- 249203**

Notified as per : G.O. No. 1002/XV-1/19-7(14)2005 AHS-1, dated 01-11-2019
Phone / Fax No. 0135 2450083 mail id : ddah.pashulok@yahoo.com

FORMET-1 : FORMET FOR SAMPLE REVIEW

FIR/Forwarding letter no :

Date :

Received from :

Sample Receipt No. :

Test requested :

Review List : (Write Yes/ No in front of each point)

- | | |
|---|---------|
| A. PROFORMA-18 from the sample submitting authority - | Yes/ No |
| B. Copy of FIR (incase FIR been filed)/ brief detail of complaint - | Yes/ No |
| C. Sample is properly sealed with lacquer - | Yes/ No |
| D. Sample not preserved/ contaminated with Formalin- | Yes/ No |
| E. Sample is properly preserved and not putrefied - | Yes/ No |
| F. Sample is in sufficient quantity - | Yes/ No |
| G. Attested copy of the Identity card of Police Personnel - | Yes/ No |
| H. Sample Seal attested by Jurisdictional Veterinary Officer - | Yes/ No |

Remarks (If any): On the basis of above report, sample is accepted/ rejected.

**Signature/ Seal of
Project Director**

Declaration by Submitting Officer:

I..... posted as
at....., hereby declare that the sample mentioned
above was opened and reviewed in my presence.

Date :

Place :

Signature

For Official Use

Sample ID generated :

FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE
SHEEP AND WOOL EXTENSION INSTITUTE
Pashulok, Rishikesh- 249203

Notified as per : G.O. No. 1002/XV-1/19-7(14)2005 AHS-1, dated 01-11-2019
Phone / Fax No. 0135 2450083 mail id : ddah.pashulok@yahoo.com

FORMET-2 : FORMET FOR SAMPLE RECEIPT

Date :

Sample Receipt No. :

Sample Description :

No. of samples :

Sample quantity and condition :

Date of sample collection :

Location of sample collection :

Name & Address of Veterinary
Officer who did the Sampling :

Name & Address of Controlling
Officer of Police Thana or
Police Chauki :

Identification details of sample
accompanying police personnel :

Test requested for :

Sample review status :

Project Director
Seal

**FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE
SHEEP AND WOOL EXTENSION INSTITUTE
Pashulok, Rishikesh- 249203**

Notified as per : G.O. No. 1002/XV-1/19-7(14)2005 AHS-1, dated 01-11-2019
Phone / Fax No. 0135 2450083 mail id : ddah.pashulok@yahoo.com

FORMET-3 : FORMET FOR SAMPLE REJECTION

Letter No.

Dated

To,

In-charge,
Police Thana/ Police Chauki
District

Sir,

Kindly refer to your FIR No. Forwarding Letter No.
dated The sample was received on dated.....
Receipt No.

The sample received with above letter is not admissible to the laboratory
because :-

1. Missing PROFORMA-18 from the sample submitting authority.
2. Missing attested copy of the Identity Card of the Sample
accompanying police personnel, attached with the forwarding letter.
3. Missing copy of FIR (incase FIR been filed) / brief detail of complaint
4. Sample has not been received in proper seal with lacquer.
5. Sample Seal attested by Jurisdictional Veterinary Officer is missing.
6. Sample has not been properly preserved.
7. Sample is too much putrefied.
8. Sample was found contaminated with Formalin.
9. Sample has not been found in sufficient quantity
10. Some other reasons narrated as

**Signature/ Seal
Project Director**

cc : 1. Veterinary Officer, Govt. Vety. Hosp. District
2. Chief Veterinary Officer, District
3. Secretary, Uttarakhand Animal Welfare Board, Top Floor, Pashudhan
Bhawan, Mothronwala, Dehradun.

**Signature/ Seal
Project Director**

FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE

FORMET-4 : FORMAT FOR SAMPLE RECEIVING REGISTER

S.No.			
Date of Receipt			
Receipt No.			
Sample quantity and condition			
Date of sample collection			
Name & Address of Veterinary Officer who did the Sampling			
Name & Address of Controlling Officer of Police Thana or Police Chauki			
Identification details of sample accompanying police personnel			
Location of sample collection			
Brief Description			
Sample ID No.			
Signature of Project Director			

**FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE
SHEEP AND WOOL EXTENSION INSTITUTE
Pashulok, Rishikesh- 249203**

Notified as per : G.O. No. 1002/XV-1/19-7(14)2005 AHS-1, dated 01-11-2019
Phone / Fax No. 0135 2450083 mail id : ddah.pashulok@yahoo.com

FORMET-5 :

FORMAT FOR SAMPLE FORWARDING to BOVINE TISSUE FORENSIC ANALYSIS LAB

Date :

Sample ID :

Quantity (approx) :

Sample preservation condition :

Test Required

- (a) Species Specific PCR (mitochondrial D Loop)
- (b) PCR by Universal Primer Technique (12s rRNA) & Species Identification by RFLP

**Signature/ Seal
Project Director**

FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE

FORMET-6 : SAMPLE ANALYSIS REGISTER OF THE LAB

			S. No.
			Date of Receipt
			Sample ID No.
			Sample quantity and condition
			Sample Description
			Test Results for Species Specific PCR (mitochondrial D Loop)
			Test Results for PCR by Universal Primer Technique (12s rRNA) & Species Identification by RFLP
			Signature of Veterinary Officer Grade-2
			Signature of Veterinary Officer Grade-1
			Signature of Project Director

**FORENSIC ANALYSIS LABORATORY FOR DETECTION OF BOVINE TISSUE
SHEEP AND WOOL EXTENSION INSTITUTE
Pashulok, Rishikesh- 249203**

Phone / Fax No. 0135 2450083 mail id : ddah.pashulok@yahoo.com

* Notified as per : G.O. No. 1002/XV-1/19-7(14)2005 AHS-1, dated 01-11-2019

FORMET-7 : TEST REPORT

Test Report No : Date :
Sample Receipt No :
Sample Receipt date :
Sample ID No :
Sample Collected by :

Forwarding Police Officer :

Brief Details about Sample :

Number of Samples :
Test Result :

Note:

- No part of report, except in full, shall be reproduced without written consent of the laboratory.
- The sample, if available, will be stored up to 1 year from the date of receipt in the laboratory unless otherwise specified.

**Signature/ Seal
Project Director**

cc : 1. Veterinary Officer, Govt. Vety. Hosp. District
2. Chief Veterinary Officer, District
3. Secretary, Uttarakhand Animal Welfare Board Dehradun.

**Signature/ Seal
Project Director**

* The lab has been Notified under Rule-33 of Uttarakhand State Protection of Cow Progeny (Amendment) Rules, 2018, read with section 30(d) of Indian Veterinary Council Act, 1984, section-45 of Indian Evidence Act, 1872 & section-293 of CrPC.